






Kartell

The automatic pipette washer is composed of

<p>Basket COD. 219 H 645 mm COD. 222 H 495 mm</p>	
<p>Jar COD. 218 H 650 mm COD. 221 H 503 mm</p>	
<p>Automatic pipette washer COD. 215 H 980 mm COD. 217 H. 734 mm</p>	
<p>Pipe attachment (internal bag with 2 pcs)</p>	
<p>Outlet tube at the base</p>	

WASHING PROCEDURE USING THE JAR

1. Place the pipette into the basket
2. Fill in the jar with the washing liquid (detergent/distilled water)
3. Leave the basket in position for sufficient time to clean the pipettes (two/three hours)
4. Remove the basket from the jar
5. Place the basket into the automatic pipette washer
6. Connect the pipe attachment provided to the automatic pipette washer (siphon connection)
7. Connect the water tap to the pipe attachment by means of a rubber or PVC tube (internal Ø 13mm not supplied)
8. Connect the outlet tube from the base to the drain

WASHING AND RINSING PROCEDURE DIRECTLY INTO THE AUTOMATIC PIPETTE WASHER

1. Fill the pipette washer with water
2. Add the detergent
3. Turn the water tap on and allow the system running until the rinsing is complete

WASHING PROCEDURE BY DISTILLED WATER DIRECTLY INTO THE AUTOMATIC PIPETTE WASHER

1. Fill the pipette washer with distilled water
2. Place the basket containing the pipettes already washed with ordinary water inside the automatic pipette washer and proceed with the rinsing by distilled water

WARNING

In case of use of chromic mixture NEVER prepare it directly in the container. Polyethylene resists up to a temperature of 80/90 ° C. It is therefore advisable to pour the cold ready made mixture, in and place the container in a 15-20 liter Polyethylene basin to avoid accidental spills of the chromic mixture.