



Dynalab® Lab

Dynalab Corp.
175 Humboldt Street, Suite 300
Rochester, New York 14610 USA

Tel 585.334.2064 800.334.7585
Fax 585.334.0241
www.Dynalab.com

Magnetic Stirrer, Non-Electric Instructions

Dynalab Part Number: 204015

1. Connect one end of the inlet hose to top connector.
2. Connect bottom outlet hose. (If water is used, run tubing to disposal source such as a sink.)
3. Connect inlet hose to pressure source (water or air). Fitting supplied for pressure source.
4. Adjust pressure as needed.

