



Disassembly/Reassembly of the Automatic Titrating Burette

Dynalon Items: 896955-0005, -0010, -0015, -0025, -0050

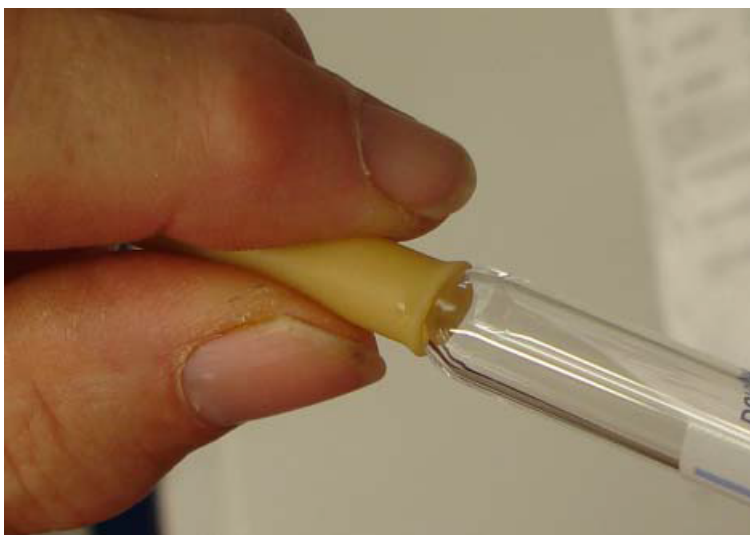


1) Remove the plastic tip from the latex hose.

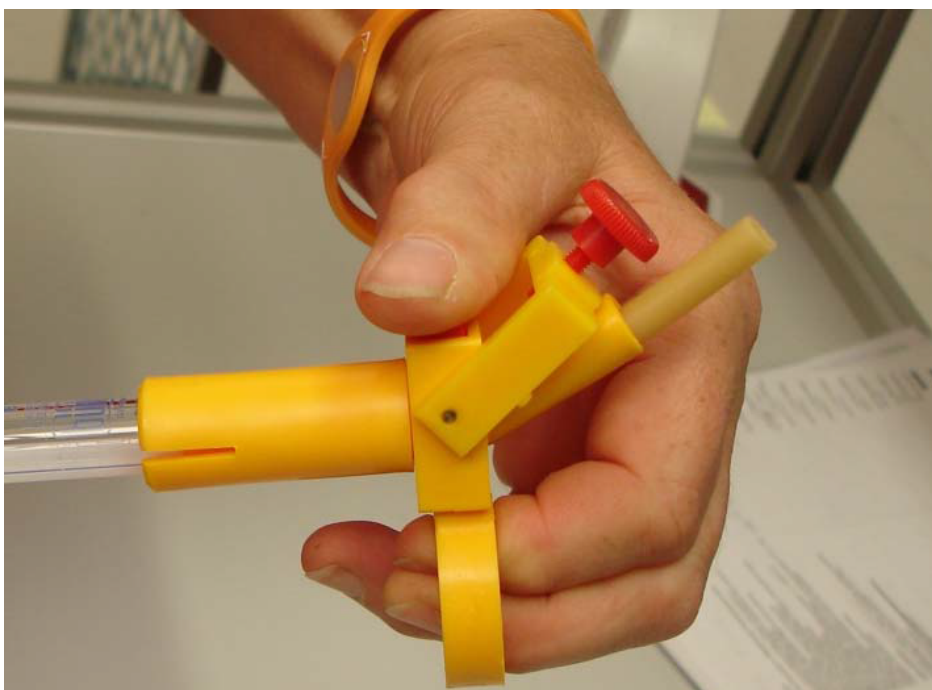

2) Push the shackle down and depress the push button all the way to take the spring pressure off the latex hose.

3) While continuing to depress the push button, slowly twist/turn and ease the burette out of the holder. The hose will come with it.

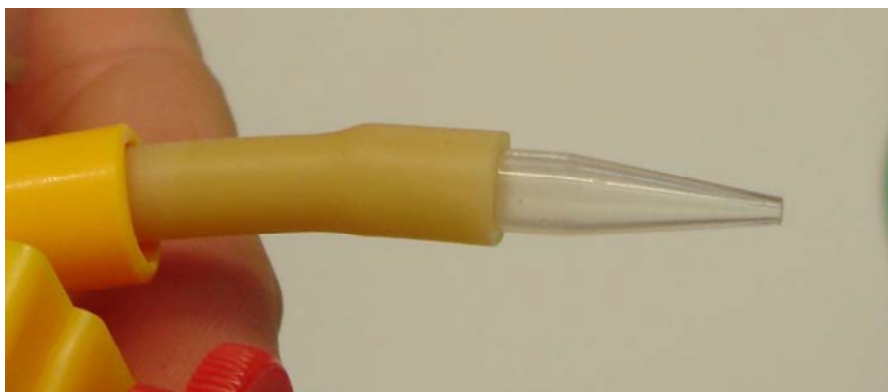

4) To reassemble, do the same steps in the opposite way.

Note: it may be useful to warm the latex or the PE tubing to remove it or to insert it onto the burette and to reinsert the plastic tip.

Cleaning:

Also, to clean the burette, it is easiest to fill the bottle with a cleaning solution and then fill the burette and let it soak. Then flush it with distilled water etc.